Safe Mixing

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Page 2

Magnetic Stirrer MR Hei-Mix S
Operating manual must be read before initial start-up.
Please follow the safety instructions provided.
Please keep for future reference.

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About this Document	4
Version and variants	4
About this manual	4
Symbols and keywords.....	5
Safety Instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / electrical safety	7
Qualifications of employees.....	7
Obligations of the operator	7
Installation site	7
Modifications to the device.....	8
Employee safety	8
California Residents	8
Safety during use	9
Disposal	9
Device Description	10
Device overview	10
MR Hei-Mix S overall view	10
Start-Up	11
Setting up device	11
Switch unit on and off	11
Operation	12
Place vessels.....	12
Start mixing	12
Set speed and start rotation.....	12

Cleaning and Maintenance	13
Cleaning.....	13
Maintenance	13
Troubleshooting	13
Possible errors and how to resolve	13
Assembly	14
Electrical connections.....	14
Disassembly and Storage	15
Dismantling, transportation and storage	15
Dismantling	15
Transportation and storage	15
Accessories and Spare Parts	15
Scope of delivery.....	15
Attachments	16
Technical data	16
China RoHS Declaration of Conformity	17
Service	19
Contact / Technical Service	19
Warranty	20
Confirmation of condition.....	21

Copyright

This operating manual is protected by copyright and is exclusively intended for the internal use of the purchaser.

Handing over this operating manual to third parties, reproducing it in any type or form – in part or in full – and utilizing and/or communicating its contents is prohibited without prior written approval from the manufacturer.

Any violations will result in an obligation to provide compensation for damages. Other claims remain unaffected.

➤ Version and variants

Version

This manual describes the function, operation and maintenance of the magnetic stirrer MR Hei-MIX S.

Version	Alteration date
1.1	10/2019

Variants

The device is available in different variants. Certain features or functions are only available in specific product variants. These variants are described in this operating manual.

➤ About this manual

This operating manual is an integral part of the device described here.

- Carefully read through this operating manual and observe all of the safety instructions and warnings notices.
- Ensure that every user has carefully read the operating manual before using the device for the first time.
- Store this operating manual so that it is accessible to all users at all times.
- Pass the operating manual on to any subsequent owner.

The current version of this operating manual can also be found at:
[https://heidolph-instruments.com/en/service/downloads/operation-manuals#Magnetic Stirrers \(MR Hei-MIX S\)](https://heidolph-instruments.com/en/service/downloads/operation-manuals#Magnetic%20Stirrers%20(MR%20Hei-MIX%20S))

In addition, also observe the local regulations at the installation site.

➤ Symbols and keywords

Standardized symbols and signal words are used in this operating manual to warn against any dangers and provide important instructions. These instructions must be strictly observed to avoid accidents and damage.

The following terms and basic symbols are used:

Symbol	Additional signal words / explanation
Warning signs 	The yellow triangle indicates hazardous situations. It is used in combination with the following signal words: DANGER: Indicates a hazardous situation which, if not avoided, will result in serious injury or death. WARNING: Indicates a hazardous situation which, if not avoided, may result in serious injury or death. CAUTION: Indicates a hazardous situation which, if not avoided, may result in property damage and minor or moderate injury.
Prohibitory signs 	Prohibited: The red circle indicates a situation that should be avoided under all circumstances and which, if not avoided, may result in serious injury or death.
Mandatory signs 	Must be observed: The blue circle indicates important information. Please observe this information to avoid any property damage.

Other symbols used:

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪	List of information
▪	
▪	
a.	List of variants
b.	
c.	

➤ General safety instructions

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

*Also see the attachment "EU Declaration of Conformity".

The device has been manufactured according to state-of-the-art technology and in compliance with recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure that the operating manual is available at all times.

The device may only be used under the following conditions:

- Only operate the device if it is in a technically perfect condition.
- Only operate the device if you are properly aware of the risks and required safety measures.
- Only operate the device in accordance with the instructions given in this operating manual.
- If any information is ambiguous or missing, ask your superior or contact the manufacturer.
- Do not operate anything on the device without authorization.
- Only use the device in accordance with its intended use.

Intended use

The device may only be operated by authorized personnel.

The device is suitable for the following uses:	<ul style="list-style-type: none"> • mixing • stirring • titrating
--	---

The device may be operated in research laboratories, other laboratories and production facilities in the following sectors:	<ul style="list-style-type: none"> • Chemistry • Pharmacy • Biology • Environmental analytics • Basic research • Similar research laboratories
---	--

Unintended use/Reasonable foreseeable misuse

Any use which deviates from the device's intended use is considered to be improper. The manufacturer is not liable for damage that occurs as a result. The risk is borne by the operator alone.

When used in medical and food sectors the operator alone is responsible for compliance with the applicable regulations and standards.

Installation / electrical safety

- The device may only be connected up if the available power supply voltage corresponds to the information stated on the rating plate for the device.
- The power supply connection must be easily accessible at all times.
- An RCD circuit breaker (residual current circuit breaker) must be fitted to the electrical system in the building so that the device is disconnected from the power circuit in the event of a fault.
- Repairs may only be performed by electricians authorised by Heidolph Instruments.
- Never operate the device with a damaged power cord.
- Always turn the device OFF and disconnect the power cord before carrying out any maintenance or repair work.

Qualifications of employees

- The device may only be operated by trained personnel.
- The device may only be operated by persons who have been instructed and supervised in its proper use by trained specialist personnel.
- The device may only be operated by specialist personnel who are above the legal minimum age.
- Other persons may only work on the device under the constant supervision of experienced and trained specialist personnel.
- This operating manual must be read and understood by all persons working with the device.
- Personnel must receive safety training that ensures responsible and safe working practices.

➤ Obligations of the operator

Installation site

- The device must be installed in a suitable location.
- The device must be installed on a firm and stable surface.
- Ensure that the device and all of its components are easily accessible at all times.
- Maintain a sufficient safety clearance to, in particular, any moving and/or hot equipment components.
- It is not permitted to place or store any objects such as accessories, tools or chemicals within this clearance area during operation.
- All screw connections must be securely tightened.
- It is not permitted to operate the device near to highly flammable or explosive substances.
- Operate the device in conjunction with an extractor hood, if working with potential harmful media (see DIN EN 14175 and DIN 12924).

Safety Instructions

- The device is designed for indoor use only and under the following ambient conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity 32 - 40 °C decreasing linearly to a maximum 50 % relative humidity
Installation altitude	0 - 2,000 m above sea level
Contamination level	2
Overvoltage category	II
Permissible supply deviations	± 10 %
- If the device is operated in corrosive atmospheres, the service life of the device will decrease based on the concentration, duration and frequency of the exposure to the corrosive atmosphere e.g. concentrated hydrochloric acid (HCl).

Modifications to the device

- It is not permitted to make any unauthorised modifications or changes to the device.
- Do not attach or install any parts that have not been approved by the manufacturer.
- Unauthorised modifications or changes will void the EC Declaration of Conformity for the device and operation of the device will no longer be permitted.
- The manufacturer is not liable for any damage, dangers or injuries that result from unauthorised modifications and changes or due to the non-observance of the instructions in this manual.

Employee safety

- Ensure that the device is only operated by qualified specialist personnel and trained employees.
- Observe the following instructions to avoid any personal injuries and property damage:
 - Laboratory regulations
 - Accident prevention regulations
 - Hazardous Substance Act
 - Other generally accepted occupational health and safety regulations
 - Local regulations

California Residents

- Important information for California residents regarding Prop 65. Please visit www.P65Warnings.ca.gov for more information.

Safety Instructions

➤ Safety during use

- Beware of the effect of magnetic fields on cardiac pacemakers or magnetic sensitive data media. Observe the according warning in chapter "Start-up", "Setting up device".
- Wear the appropriate protective clothing when working on the device (protective glasses and, if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not carry out work with naked flames in the vicinity of the device (risk of explosion).
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range ($3 \cdot 10^{11}$ Hz to $3 \cdot 10^{15}$ Hz).
- Do not operate or assemble appliances in the vicinity of the device which constitute emission or radiation sources for ionizing radiation or in the ultrasonic range.
- Do not operate the device where adiabatic compression or shock waves might occur (shock wave combustion).
- Do not use substances, where the energy input due to mixing might pose a hazard.
- Do not spill liquids over the device or any parts of it.
- Remove any accidentally spilled liquids immediately.
- Eliminate errors immediately.
- Always switch the device OFF after use.
- Do not use abrasive material to clean the surface. Only wipe with damp cloth.

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device in accordance with the relevant national regulations.
- Dispose of the packaging material in accordance with the appropriate national regulations.

➤ Device overview

MR Hei-Mix S overall view

➤ Setting up device

Warning:
Magnetic field!

The magnetic field beneath the top plate could cause irritations on your cardiac pacemaker, if you get too close.
Personnel with a cardiac pacemaker should keep a minimum distance of 10 cm (4 inches) to the device.

Warning:
Risk of slipping device!

As a result of vibrations generated during operation the device might slide off the table top.
The surface must be smooth, clean and temperature resistant and should be properly leveled.

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

➤ Switch unit on and off

The ON/OFF switch is situated on the front panel.

Switch unit on

- Press the ON/OFF switch on the lower side.
- ✓ The unit is switched on.

Switch unit off

- Press the ON/OFF switch on the upper side.
- ✓ The unit powers off.

ON/OFF switch off

➤ Place vessels

Caution:
Risk of breakage and spillage

If the vessels are not placed securely they may slip off and spill the sample. Prior to start mixing ensure that all vessels are securely placed.

Accessories for round flasks and beakers are available.

- To place a number of vessels securely on plate use optional adaptors (see general catalogue).
- Single vessels with flat bottoms like beakers may be placed directly on magnetic plate.

A single flask should be arranged in the middle of the shaker plate. Several flasks should be distributed equally on the plate.

➤ Start mixing

Warning:
Risk of injury!

Returning power after a power cut off will restart mixing mode. There is a risk of injury in the vicinity of moving parts of device.

Before working with the device or in the vicinity ensure all controls are switched to "0" or "OFF" mode, when device is passiv.

When using device inside conditioning cabinets, make reference to ambient conditions as stipulated in chapter "Technical data".

Set speed and start rotation

Warning:
Risk of poisoning!

Open vessels and too high speed may result in samples splashing.

Use vessels with narrow neck especially if sample is dangerous or toxic.

Adjust speed step by step until you have reached the required rpm settings.

Wear safety glasses and especially with dangerous and toxic samples appropriate protective clothing.

Place vessel with sample and stirring bar on magnetic plate.

- Turn item on with ON/OFF switch.
- Set speed with analog knob.
- ✓ Rotation starts immediately.

➤ Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

Warning:
Risk of electric shock!

If the device is still connected to power during cleaning, you might get in contact with live parts due to accidentally entering liquids.

Prior to cleaning:

- ensure that the device is switched OFF
- ensure that the plug is pulled from the electric socket.

Caution:
Damage to the device

Accidentally entering liquids will cause the device to fail .
Only wipe with a damp cloth. Never use a soaked cloth.

Attention:
Risk of damage to the surfaces

All of the surfaces may be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent containing chlorine
- Solvent-bearing agents (e.g. acetone)
- Ammonia
- Abrasive cleaning agents such as scouring agents, cleaning wool or agents with metallic particles

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

➤ Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

➤ Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
No rotation function	No power	<ul style="list-style-type: none"> ▪ Check size and shape of plug and compatibility to your electrical socket ▪ Check circuit breakers
	No magnetic bars in vessel rpm set to "0"	Insert magnetic bars Set speed

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

Electrical connections

Repairs of any kind may only be carried out by a **qualified electrician certified by Heidolph Instruments**. Improper repairs may result in considerable danger. Contact your local Heidolph Instruments distributor if repairs are required.

Warning:
Electric shock!

The unit and voltage must match. The type label on the back side of the unit provides all voltage specifics.

Rating plate

Color code for electrical connections:

Europe

BLUE

N: Neutral conductor

BROWN

P: Phase

North America

WHITE

N: Neutral conductor

BLACK

P: Phase

The device is connected to the mains with the power cord supplied. This plugs into the socket on the rear of the housing.

For countries where a different plug than the standard plug supplied is used:

- The plug supplied may only be changed by a professional electrician.
- If you wish to use an adapter, ensure that it is grounded and approved in accordance with the local regulations.
- Wiring, adapter and plug must at least meet the electrical power stated on the rating plate.

➤ Dismantling, transportation and storage

Dismantling

**Warning:
Risk of injury!**

While the unit is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling unit ensure that the unit is switched off and the plug is pulled from the electric socket.

- ✓ The unit is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the unit.
- Remove all vessels.
- Disassemble attachments.

Transportation and storage

- Store and transport the unit and its components only if they were emptied, cleaned and dismantled.
- To carry the device wear safety gloves and lift the unit from below.
- Store and transport the device and parts of device in the original packing material or alternatively in an adequate container to prevent damage.
- Seal the package carefully against unauthorized or accidental opening.
- Store the device in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Quantity	Product number
MR Hei-Mix S*	1	503-02000-00
Operating Manual English / German	1	01-005-004-60
Warranty registration / Confirmation of condition	1	01-006-002-78
EC Declaration of Conformity	1	01-001-025-02
Power cord (fixed)	1	depending on country

* The product number is valid for 230 V devices in the european community. Please contact Heidolph Instruments for further product numbers on country specific varieties.

➤ Technical data

MR Hei-Mix S

Standard supply voltage	230 V/50 Hz or 115 V/60 Hz
Supply power (W)	7
Protective class (IEC 61140)	2 <input type="checkbox"/>
Protection class (IEC 60529)	IP 21
Drive	Shaded pole motor
Operating mode	continuous
Speed range (rpm)	0 - 2,200
Speed accuracy (%)	±5
Stirring capacity, max (H ₂ O) (l)	5
Load capacity, max. (kg)	6
Plate diameter (Ø) (mm)	104
Plate material	PVDF
Temperature resistance max (°C)	40 (104 °F)
Weight (kg)	1.1
Dimensions (l x w x h) (mm)	140 x 126 x 80

China RoHS DECLARATION OF CONFORMITY

Heidolph Instruments GmbH & Co.KG has made reasonable efforts to ensure that hazardous materials and substances may not be used in its products.

In order to determine the concentration of hazardous substances in all homogeneous materials of the subassemblies, a "Product Conformity Assessment" (PCA) procedure was performed. As defined in GB/T 26572 the "Maximum Concentration Value" limits (MCV) apply to these restricted substances:

- Lead (Pb): 0.1%
- Mercury (Hg): 0.1%
- Cadmium (Cd): 0.01%
- Hexavalent chromium (Cr(VI)): 0.1%
- Polybrominated biphenyls (PBB): 0.1%
- Polybrominated diphenyl ether (PBDE): 0.1%

Environmental Friendly Use Period (EFUP)

EFUP defines the period in years during which the hazardous substances contained in electrical and electronic products will not leak or mutate under normal operating conditions. During normal use by the user such electrical and electronic products will not result in serious environmental pollution, cause serious bodily injury or damage to the user's assets.

The Environmental Friendly Use Period for Heidolph Instruments GmbH & Co.KG products is 25 years.

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This table is created according to SJ/T 11364-2014

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Note: Table applies to all products. Some of the components or parts listed above may not be part of the enclosed product.

- O: 表示该有毒有害物质在该部件所有均质材料中的含量均在GB/T 26572规定的限量要求以下。
- O: Indicates that the above mentioned hazardous substance contained in all homogeneous materials of the part is below the required limit as defined in GB/T 26572.
- X: 表示该有毒有害物质至少在该部件某一均质材料中的含量超出GB/T 26572规定的限量要求。
- X: Indicates that the above mentioned hazardous substance contained in at least one of the homogeneous materials of this part is above the required limit as defined in GB/T 26572.

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Apart from the disclosures in the above table, the subassemblies are not intentionally manufactured or formulated with lead (Pb), mercury (Hg), cadmium (Cd), hexavalent chromium (CrVI), polybrominated biphenyls (PBB), and polybrominated diphenyl ethers (PBDE).

Products manufactured by Heidolph Instruments GmbH & Co.KG may enter into further devices or can be used together with other appliances.

With these products and appliances in particular, Heidolph Instruments GmbH & Co.KG will not take responsibility for the EFUP of those products and appliances.

Place, date Schwabach, 19.09. 2019

Wolfgang Jaenicke
Chief Executive Officer CEO

Marcell Sarré
Vice President
Quality Management &
Technical Service

➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning! **Danger of poisoning!**

Contaminated devices can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
- take adequate protective action towards our parts receiving and service personnel
- mark the pack in accordance with Ordinance on Hazardous Substances

A "Confirmation of Condition" form can be found at the end of this operating manual.

→ Prior to shipping the device for repair, complete a copy of this form and submit it in advance.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 – 9122 - 9920-74
Fax: +49 – 9122 - 9920-84

E-Mail: service@heidolph.de
www.heidolph.com

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at
www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3 year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty period begins with the date of purchase. The serial number is also valid without registration.

This warranty covers material and manufacturing defects.

In the event of a material or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damage incurred as a result of improper handling or transport.

Warranty claim?

- ➔ Please notify Heidolph Instruments or your local Heidolph distributor if you wish to make a warranty claim.

➤ Confirmation of condition

→ In the case of repair, copy and complete the Confirmation of condition form and send it to Heidolph Instruments.

1. Details about the device

Product number _____

Serial number _____

Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the device in a condition which does not represent any health threats for the staff of our service department?

Yes _____ No _____

If not, which substances has the device come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date

Signature

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG